

Technical Data Sheet

Eye Wash Unit Aqua Push TWIN



Order no.: 177 525
EAN: 4013207177525

Product designation:

Eye Wash Unit Aqua Push TWIN, with 2 shower heads

The Eye Wash Unit Aqua Push TWIN is manufactured according to factory specifications and complies with DIN EN 15154-2 and the BGI/GUV-I 850-0 "Safe working in laboratories - principles and guidelines" issued by the German Social Accident Insurance (DGUV).


Eye showers with a water connection are mandatory first aid equipment for workplaces where hazardous substances are handled. The eye shower is used for targeted rinsing of the eyes and for rinsing the face



DESCRIPTION

- Practical emergency eyewash featuring two pearl-shaped spray heads for simultaneous flushing of both eyes as an immediate first-aid measure
- Ergonomically designed plastic handle with an integrated locking mechanism for easier, hands-free operation during rinsing
- The specially designed soft spray heads of the Aqua Push TWIN Eyewash provide an effective, wide-coverage water spray
- Adjustable water flow directly on the eyewash unit by continuously turning the flow control valve to suit individual requirements
- At the fully open flow setting, the flow rate is approximately 7 L/min per spray head
- Automatically opening dust covers protect the eyewash unit from contamination when not in use
- Supplied with a high-quality 150 cm stainless steel braided hose and a standard ½ inch (ISO 228) connection
- Quick and easy installation by simply connecting the unit to the local cold drinking water supply
- Included also: a mounting bracket for convenient tabletop installation and a self-adhesive safety sign for standards-compliant identification

MAIN FEATURES

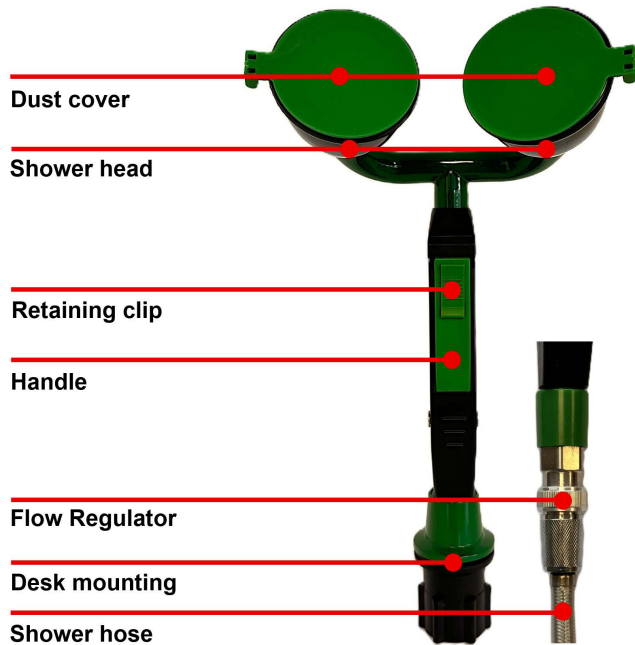

MATERIALS

The materials used are particularly kind to the skin and therefore have no negative effect on human health

- | | |
|-----------------|---------------------------------------|
| • Handle: | Acrylonitrile-Butadiene-Styrene (ABS) |
| • Shower head: | High-density polyethylene (HDPE) |
| • Dust cover: | Acrylonitrile-Butadiene-Styrene (ABS) |
| • Deskmounting: | Acrylonitrile-Butadiene-Styrene (ABS) |
| • Shower hose: | Stainless steel braided hose |

WEIGHT

- approx. 1020 g

MARKING

- Trade mark of manufacturer (EKASTU Safety) on left dust cover
- Order number and Made in Germany on right dust cover

DONNING AND USING

- Store the Eye Wash Unit Aqua Push TWIN and spare parts in a dry place, protected from dust and corrosive and oxidant agents

HANDLING

- The eye shower should be connected to a cold drinking water connection.
- The location of the eye shower must be known to all employees
- Please consider the information brochure!

USER REFERENCES

EKASTU Safety GmbH guarantees the indicated achievement according to class and type. It is to be noted that the laboratory test values can considerably deviate from those, which are reached in practice. The user must read and understand all functional information. Use the EKASTU-Eye Wash Unit Aqua Push with this information brochure, the relevant valid statutory regulations and the safety requirements of the profession associations for example DGUV Information 204-022 "First Aid in Operation".

TD 177 525 English_01-07.26